# BAL series

ES | EN | FR

V.1.2  
05/08/2024

BALANZA ANALÍTICA VERIFICADA

VERIFIED ANALYTICAL BLANCE

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# EN

## INDEX

<b>1. INSTALLATION INSTRUCTION</b>	<b>45</b>
<b>2. STORAGE</b>	<b>46</b>
<b>3. WEIGHING PAN ASSEMBLY</b>	<b>46</b>
<b>4. KEYBOARD AND DISPLAY</b>	<b>47</b>
<b>5. OVERVIEW</b>	<b>47</b>
5.1 FRONT PART OF THE BALANCE, M/MW/ L / LW MODELS	47
5.2 REAR AND BOTTOM PART OF THE BALANCE	48
5.3 REAR AND BOTTOM PART OF THE BALANCE	49
<b>6. WEIGHING</b>	<b>50</b>
6.1 STAND BY	50
6.2 SIMPLE WEIGHING	50
<b>7. CALIBRATION</b>	<b>50</b>
7.1 EXTERNAL CALIBRATION	50
7.2 INTERNAL CALIBRATION	51
7.2.1 AUTOCALIBRATION (AUT-CAL)	52
7.2.2 INTERNAL CALIBRATION (I-CAL) [LOCKED FOR VERIFIED BALANCES]	52
7.2.3 EXTERNAL CALIBRATION (E-CAL) [LOCKED FOR VERIFIED BALANCES]	52
7.2.4 TECHNICAL CALIBRATION (TEC-CAL) [LOCKED FOR VERIFIED BALANCES]	52
7.3 VERIFICATION	53
<b>8. TARE FUNCTION</b>	<b>54</b>
8.1 MANUAL TARE FUNCTION	54
<b>9. WEIGHT UNITS</b>	<b>55</b>
<b>10. PC COMMUNICATION SETTING</b>	<b>56</b>
<b>11. PRINTER SERIAL COMMUNICATION SELECTION</b>	<b>57</b>
<b>12. TRANSMISSION SPEED SELECTION</b>	<b>57</b>
<b>13. AUTOZERO FUNCTION</b>	<b>58</b>
<b>14. FILTERS SELECTION</b>	<b>59</b>
<b>15. STABILITY FUNCTION</b>	<b>59</b>
<b>16. DISPLAY CONTRAST REGULATION</b>	<b>60</b>
<b>17. DISPLAY BACKLIGHT SETUP</b>	<b>60</b>
<b>18. AUTO POWER-OFF FUNCTION</b>	<b>61</b>
<b>19. SELECTION OF WORKING MODE WITH TABLET</b>	<b>62</b>
<b>20. PIECECOUNTING FUNCTION</b>	<b>63</b>
20.1 VISUALIZATION OF TOTAL AND UNIT WEIGHT OF PIECES	64
20.2 VISUALIZATION OF TOTAL AND UNIT WEIGHT OF PIECES	64
20.3 AUTOMATIC UPDATE OF UNIT WEIGHT	65

<b>21. DENSITY DETERMINATION OF A SOLID OR A LIQUID</b>	<b>65</b>
21.1 DENSITY DETERMINATION OF A SOLID OR A LIQUID	65
21.2 DENSITY DETERMINATION OF A LIQUID	66
<b>22. MAXIMUM LOAD DETERMINATION FUNCTION</b>	<b>67</b>
<b>23. PERCENTAGE WEIGHING FUNCTION</b>	<b>68</b>
23.1 MODE WITH WEIGHT REFERENCE	68
23.2 MODE WITH MANUAL INSERTION OF THE REFERENCE WEIGHT	68
<b>24. ANIMAL WEIGHING FUNCTION</b>	<b>69</b>
<b>25. TOTALIZING</b>	<b>70</b>
<b>26. THRESHOLD FUNCTION</b>	<b>71</b>
26.1 BOTH THRESHOLDS SET	71
26.2 LOWER THRESHOLD SET ONLY	71
26.3 HIGHER THRESHOLD SET ONLY	72
<b>27. RS232 INTERFACE FEATURES</b>	<b>72</b>
<b>28. ERROR CODES</b>	<b>76</b>
<b>29. MAINTENANCE AND CARE</b>	<b>76</b>
<b>30. QUICK GUIDE TO BALANCE PARAMETERS SETUP</b>	<b>77</b>
<b>31. QUICK GUIDE TO THE USE OF THE BALANCE PROGRAMS</b>	<b>78</b>
<b>32. BALANCE TECHNICAL CHARACTERISTICS</b>	<b>79</b>
<b>33. WARRANTY</b>	<b>79</b>
<b>34. EQUIPMENT DISPOSAL</b>	<b>79</b>

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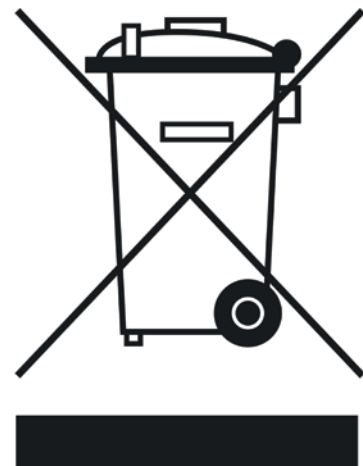
# 1. INSTALLATION INSTRUCTION

## WARNING:

Please follow carefully these steps for installing and use the new balance before starting your work routine. A way of use of the instrument different from this user manual will not guarantee the instrument's safety anymore.

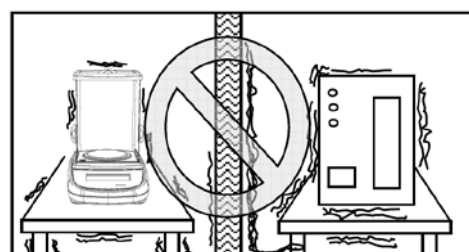
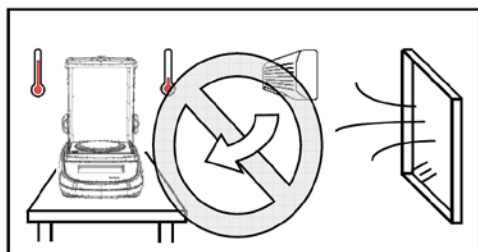
- **Remove** the balance, calibration weight( if included) and accessories from the cartoon and check for any visible damage of the instrument.



## ATTENTION

In models with internal calibration remove the screw in the left lower part of the balance (see par.5).

- **Do not install** the balance in a place with air flows, heavy thermal changes and vibrations.
- **Do not use** the balance in blast risk environment.
- **The humidity rate** of the balance environment must be between 45% and 75%..

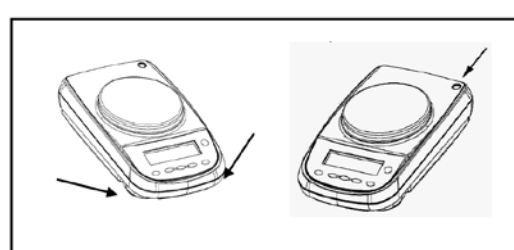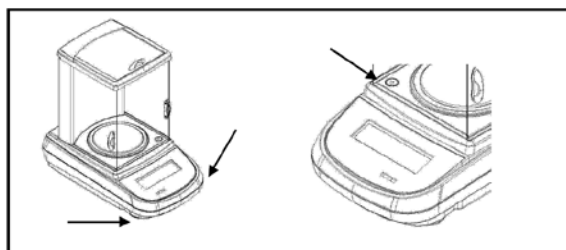- **Place** the weighing pan and the support pan on the balance (see par. 3).
- **Level** the balance using the level bubble and levelling feet located underneath the case.

- **Connect** power supply to connector 2 located on the rear panel of the unit (see par. 5).
- **Insert** power unit into AC outlet, **which shall be easily accessible**; after few seconds the balance will automatically switch on.
- Balance Warm-up:
  - Wait 8 hours from switch one for Analytical balanaces (0,0001g).
  - Wait 30 minutes for other models.

Then calibrate the balance using the appropriate mass (if supplied), following instructions (par.7).


**It is strongly suggested not to disconnect the balance from the electrical mains and to use the  button to stand-by the balance when you finished using it.**

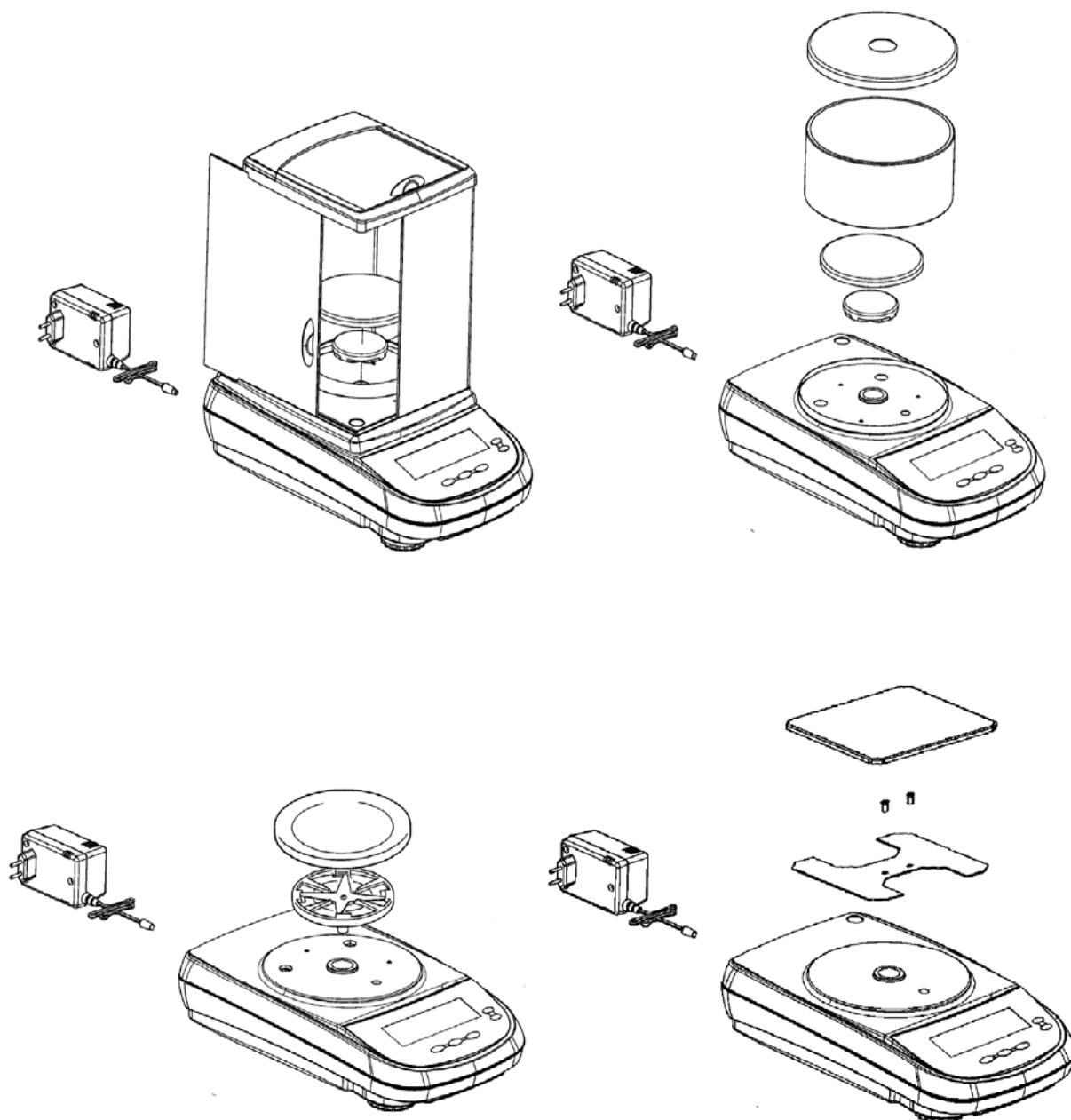- **Calibrate** again the balance every time it's moved to another place.
- **Check** balance calibration periodically.
- **We recommend** not dropping heavy objects on balance pan, in order to avoid damages.
- **Assistance service** must be effected by specialized staff and the spare parts used must be original. Therefore, it is necessary to apply to the reseller who sold the equipment.


## 2. STORAGE

- Storage temperature: +5 °C... +40°C
- Storage humidity 45% - 75%.
- Keep package in the case the balance should be sent back to the factory for service. Disconnect the cables and the accessories for avoiding damages during transportation.
- Do not place the balance in extreme temperature and humidity condition and avoid the balance to take strong hit.

## 3. WEIGHING PAN ASSEMBLY


## 4. KEYBOARD AND DISPLAY
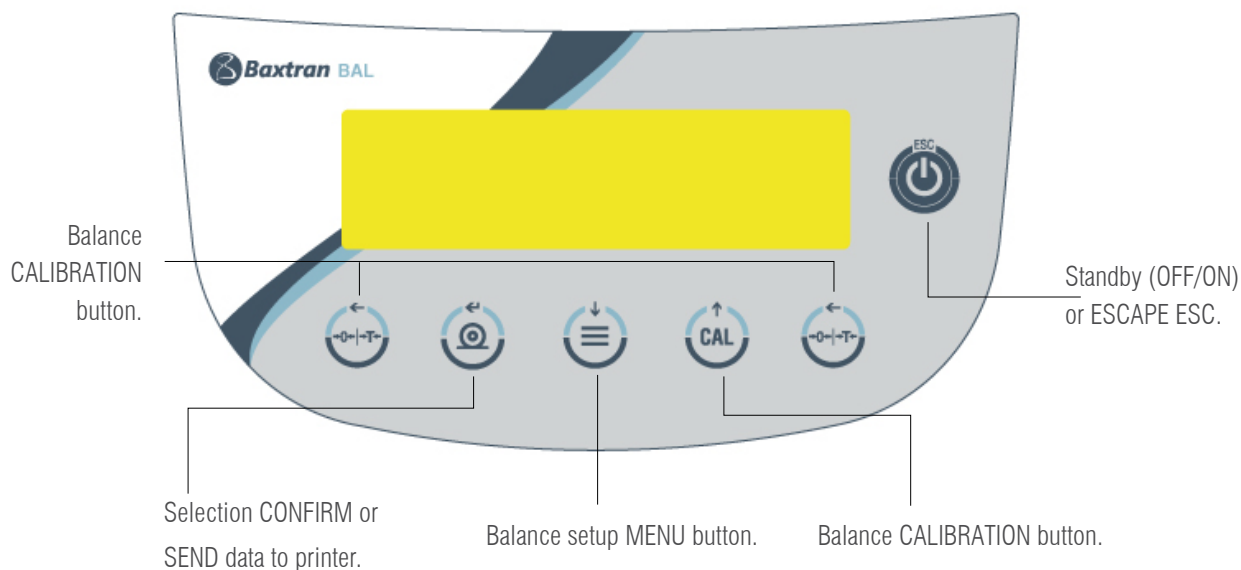
<b>*</b>	Stability indicator.	<b>▼</b>	Insert data mode.
<b>0</b>	Zero indicator.	<b>H</b>	Upper threshold.
<b>%</b>	Percentage weighing.	<b>L</b>	Lower threshold.
<b>PC</b>	Piece counting.	<b>DS</b>	Density measure mode.
	Battery charge indicator.	<b>ct, ozt, lb, GN, dwt, kg, mg</b>	Measure unit

## 5. OVERVIEW


### 5.1 FRONT PART OF THE BALANCE

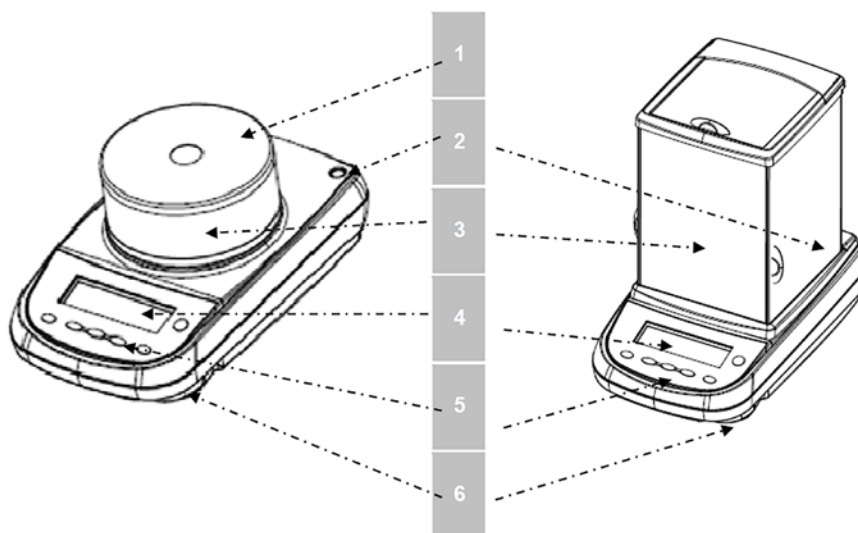**Fig.1 Balance with circular and three-panes windshield**

- |                      |                     |
|----------------------|---------------------|
| 1. Cap of windshield | 4. LCD display      |
| 2. Levelling bubble  | 5. Function buttons |
| 3. Windshield        | 6. Levelling feet   |


## 5.2 REAR AND BOTTOM PART OF THE BALANCE

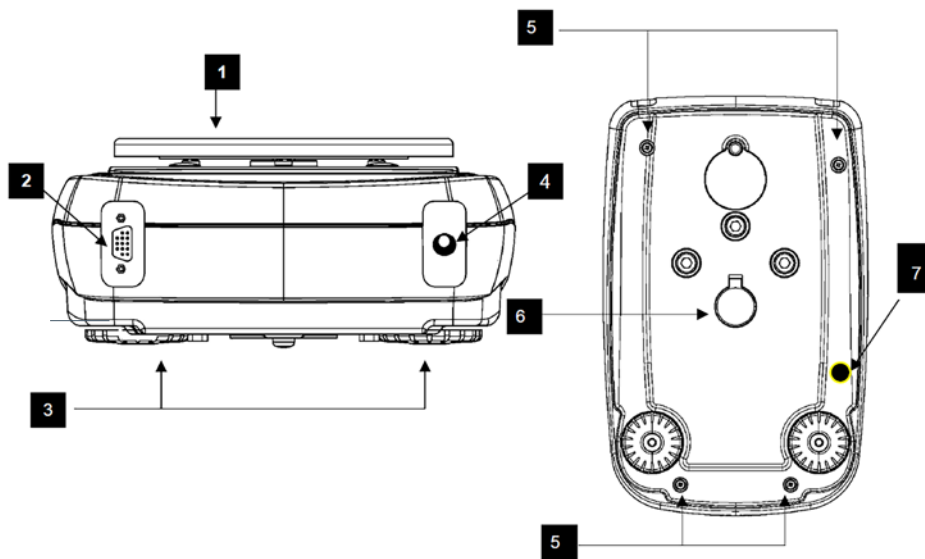**Fig.2** Rear and bottom part of the balance

1. Weighing pan	6. Hook for under-balance weighing. Hook not available in internal calibration models and for verified balances. <b>Note:</b> please remove the cap to access the hook
2. Nine poles (pin) female connector for RS232 interface for printer/PC/external keyboard	
3. Levelling feet	7. Protection screw for autocalibration motor (only for models with internal calibration). <b>ATTENTION:</b> remove this screw by hand as soon as the balance is removed from its package
4. connector for power supply	
5. Screws for balance closing. In models with 4 feet it is necessary first to remove the two rear feet (one fixed and one moving) to access to the closing screws	

### 5.3 REAR AND BOTTOM PART OF THE BALANCE SQUARE PAN VERSION

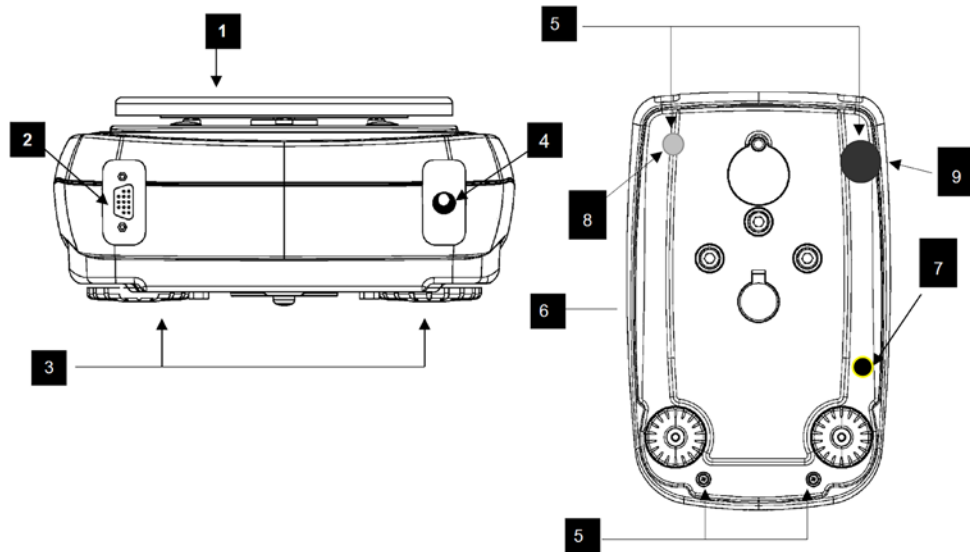**Fig.3** Rear and bottom part of the balance

1. Weighing pan	6. Hook for under-balance weighing. Hook not available in internal calibration models and for verified balances. <b>Note:</b> please remove the cap to access the hook
2. Nine poles (pin) female connector for RS232 interface for printer/PC/external keyboard	
3. Levelling feet	7. Protection screw for autocalibration motor (only for models with internal calibration). <b>ATTENTION:</b> remove this screw by hand as soon as the balance is removed from its package
4. connector for power supply	
5. Screws for balance closing. In models with 4 feet it is necessary first to remove the two rear feet (one fixed and one moving) to access to the closing screws	8. Adjustable rear
	9. Fixed rear feet

## 6. WEIGHING

- i** For verified Balances models at switch on it will appear *WARM UP* on display and countdown will start.

After having connected the balance to AC outlet, it will perform an internal circuits test, after that the balance will set itself in stand-by mode.

### 6.1 STAND BY

From "STAND BY" mode:

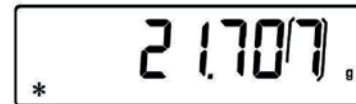
- Press  button to bring balance to work conditions.
- Press again  button to return to "STAND BY" condition.

### 6.2 SIMPLE WEIGHING

From "STAND BY" mode:

Load the sample to weigh on the pan and read the value on display as soon as the stability symbol \* (star) appears.

- i** The digit that represents the verification division "e" is rounded with a frame.

## 7. CALIBRATION

Electronic balances take mass measurements making use of gravity (g). Difference of latitude in geographic areas and altitude will vary gravity acceleration value (g). Therefore, for accurate measurements, the balance must be adjusted to the local environment. This adjustment is accomplished by calibration function.

### 7.1 EXTERNAL CALIBRATION

- i** For verified Balances models the External calibration function is locked.

Calibration is accomplished by pressing  button.

1. Press  button when pan is empty, dashes are displayed on the display.

2. When calibration weight value starts to flash, load the weight on the pan.

3. The display will stop flashing, indicating calibration weight value.

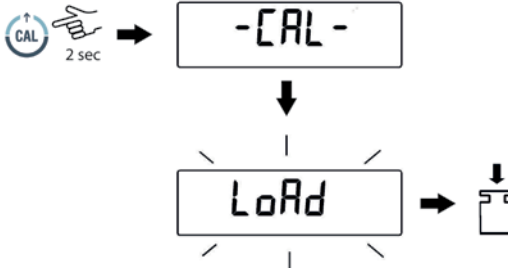
Once the calibration is effected will be shown the value of the calibrated weight and the current unit of measure.

4. Unload calibration weight from the pan.  
The balance is ready for weighing operations.

**NOTE:** if there is interference during calibration process, an error message will be displayed.

Moreover, it is possible to calibrate the balance with a calibration weight higher than the one set by default:

1. Press and keep  button pressed with empty pan until the acoustic alarm stops, then release the button. On display it will be visualized the string **-CAL-**, followed by flashing string **LoAd**

2. Load on the pan a weight equal higher or lower than default calibration weight; the balance will recognize it as valid weight if equal or higher than calibration weight as long as it is a whole number in comparison with the most meaningful digit of calibration weight .

*e.g.: if calibration weight is 200g, it will be possible to calibrate the balance with values from 100g 200g, 300g, 400g up to the highest limit of balance weighing range.*

The message **LoAd** on display will stop flashing. Once calibration has been effected, the value of calibrated weight will be displayed.

3. Unload calibration weight from the pan.  
The balance is ready for weighing operations.

**NOTE:** if there is interference during calibration process, an error message will be displayed.

## 7.2 INTERNAL CALIBRATION

**\* only for models with internal calibration “-i”**

In these balance models there are 4 calibration modes:

From display zero condition, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message **units** will be visualized on display, press then  button until you visualize **Calib** on display. Press  to confirm..

1. Select the calibration mode you wish by pressing  button in sequence:

- **AUT-CAL:** auto calibration
- **I-CAL:** internal calibration\*
- **E-CAL:** external calibration\*
- **TEC-CAL:** technical calibration\*

\*Locked for verified balances models

2. Press  button to confirm “**AUT-CAL**”, “**I-CAL**”, “**E-CAL**”.

To confirm “**TEC-CAL**” keep pressed the  button until the acoustic alarm gets mute.

3. After selection, the balance returns to calibration menu. Press and keep pressed  button until the acoustic alarm gets mute, then release the button. Balance is again ready for weighing operations

## 7.2.1 AUTOCALIBRATION (AUT-CAL)

### This is Factory setting for verified balances models

The balance self-calibrates when the temperature variation exceeds the factory preset value and at factory preset time intervals, through the internal reference mass, and only if the balance pan is empty.

When the balance needs to perform the Automatic calibration, the display will show the following message: *CAL 25*

A 25-seconds countdown will start during which you can decide if:

- Stop the automatic-calibration procedure by pressing the  button that will be delayed of 5 minutes.
- Let the countdown finish so that the automatic calibration starts.

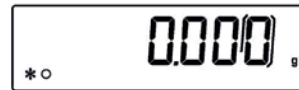
**Note:** during the countdown do NOT load anything on the plate!

In this mode, it is also possible to carry out the calibration with the internal reference mass by pressing the  button at any moment, first ensuring that no weight is loaded on the plate.

1. Press  button with empty pan. The message *--CAL--* will be displayed and balance calibration will be effected automatically.

2. At the end of calibration, balance returns to normal weighing conditions.

If calibration is not finished due to vibrations or air flows, then the message *CAL BUT* is displayed. Press again  button. If the problem repeats then select external calibration and contact the product supplier.

## 7.2.2 INTERNAL CALIBRATION (I-CAL) [LOCKED FOR VERIFIED BALANCES]

The balance calibrates itself with internal reference mass **ONLY** on user command by pressing  button. Before effecting internal calibration, be sure that there's no weight loaded on the balance pan.

## 7.2.3 EXTERNAL CALIBRATION (E-CAL) [LOCKED FOR VERIFIED BALANCES]

The balance is calibrated through an external reference mass (**NOT** provided with models with internal calibration).

(Follow procedures described at paragraph 7.1)

## 7.2.4 TECHNICAL CALIBRATION (TEC-CAL) [LOCKED FOR VERIFIED BALANCES]

This function allows storing the value of internal reference mass whenever checking or assistance actions require it.

1. After having selected the **TEC-CAL** calibration mode, press CAL button at empty pan. It will be displayed *--CAL--*

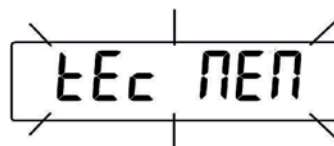
2. When the value of calibration weight start flashing on display, load the weight on to the balance pan.

3. Wait the acoustic alarm and that the displayed calibrated weight value stops flashing, then unload the weight from balance pan.

4. When string *0.000* is displayed continuously, then press and keep pressed the  button. This starts the internal weight value automatic acquisition and store. During the acquisition cycle, the display will show *tEe nEN*

5. After having stored the value of internal calibration weight, balance returns to normal weighing conditions.

6. Return to calibration menu as described at paragraph 6.2 and set the desired calibration mode: internal, automatic or external.

**ATTENTION**

**This procedure must be effected only using E2-class reference masses.**

## 7.3 VERIFICATION

General introduction:

According to EU directive 90/384/EEC balances must be verified if they are used for legally controlled area as follows:

- a) For commercial transactions if the price of goods is determined by weighing.
- b) For the production of medicines in pharmacies as well as for analyses in the medical and pharmaceutical laboratory.
- c) For official purposes.
- d) For manufacturing final packages.

In cases of doubt, please contact your local trade in standard.

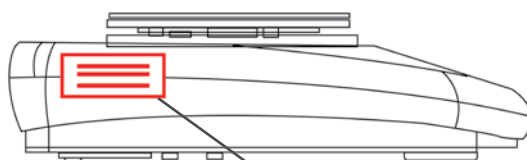
### Verification instructions

An EU type approval exists for balances described in their technical data as verifiable. If a balance is used where obligation to verify exists as described above, it must be verified and re-verified in regular intervals.

Re-verification of a balance is carried out according to the respective national regulations. The legal regulation of the country where the balance is used must be observed!

After verification the balance is sealed at the indicated positions.

**Verification of the balance is invalid without the “seal”**


Position the seal for certified balances

### **Balances with obligation to verify must be taken out of operation if:**

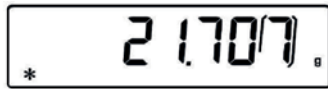- The weighing result of the balance is outside the error limit. Therefore, in regular intervals load balance with known test weight (ca. 1/3 of the max. load) and compare with displayed value.
- The re-verification deadline has been exceeded.

## 8. TARE FUNCTION

1. Load the container on the pan. The display will show the weight.


2. Press  button.  $0-t$  string will be displayed.

3. After reaching stability, the value  $0.000$  will be displayed.  
If the stability is not reached (due to air flows or vibrations or other disturbs) the dashes will remain displayed.

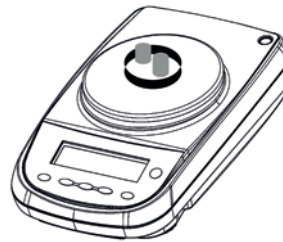4. Load the objects to weigh in the container. Read net weight value on display.

### 8.1 MANUAL TARE FUNCTION

This function allows to insert manually the tare values.

1. Press and keep pressed the  key with no objects on the weighing pan until the beep alarm stops, then release the key.

2. Will be shown on the display the following string:

3. Insert now the desired tare value using the keys  and  for increase or decrease the value, and press the  key for skip to the following digit. During the value insert mode, if keep pressed the  key is possible to delete the inserted value.

4. Press  key to confirm the value inserted.

## 9. WEIGHT UNITS

The balance can be set to display the weight in the different units, one primary (**unit1**) and one secondary (**unit2**).  
When we supply the balance, the default unit of measurement is **unit1**.

1. From display zero condition press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *unit 1* will be displayed. Then press  to confirm.
- 
2. It will be displayed *GrAN*.unit. Pressing the  or  button, it will be possible to scroll forward or backward the weight units menu.

3. Press  button to confirm or  button to shift to the other weight unit.

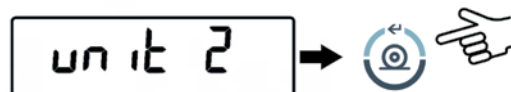
SYMBOL	UNIT	CONVERSION FACTOR 1g =
<i>GrAN</i>	GRAM	1.
<i>CARAt</i>	CARAT	5.
<i>ounCE</i>	ONCE**	0.035273962
<i>Pound</i>	POUND**	0.0022046226
<i>PEnn.</i>	PENNYWEIGHTS**	0.643014931
<i>onCEtr.</i>	ONCE TROY**	0.032150747
<i>GrANo</i>	GRAIN**	15.43235835
<i>LAEL Hon</i>	HONG KONG TAEI **	0.02671725
<i>* LAEL SGP</i>	SYNGAPORE TAEI **	0.02646063
<i>* LAEL roC</i>	R.O.C. TAEI **	0.02666666
<i>* NoNNE</i>	MOMME**	0.2667
<i>n 10</i>	x 10**	10.
<i>n 100</i>	x 100**	100.

\*\*not available in verified balances

4. After setting **unit1** (by pressing the  button to confirm), press the  button to select the second unit of measurement.

5. The "**unit2**" will be displayed, then press the  key to confirm.

6. The *GrAN* unit will be displayed. Pressing the  or  key will now be possible to scroll backwards or forwards the secondary units menu

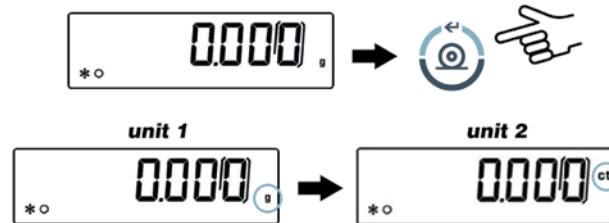
7. Press the  key to confirm or  to change to another unit of measurement (the units of measurement available are the same as those listed in point 3).

8. To escape from parameters setup menu, press the  button until the acoustic alarm gets mute, then release the button.

9. The balance returns to normal weighing conditions.

It is useful to set a second unit of measure when it is necessary to quickly display the result of a weighing in two different units

10. By setting both units of measurement, returns to normal weighing condition, it will be sufficient to press the  button until the acoustic alarm gets mute, then release the button to switch from one unit of measurement to another.

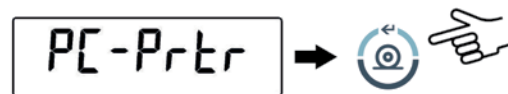
**NOTE:** Putting the balance in **Stand-by** status using the  button, the weight in the last selected measurement unit will be displayed when the power is turned on again.

Instead, by disconnecting the instrument from the electrical network, when it is switched on again, the weight will be displayed in the unit of measure corresponding to **unit1**.


## 10. PC COMMUNICATION SETTING

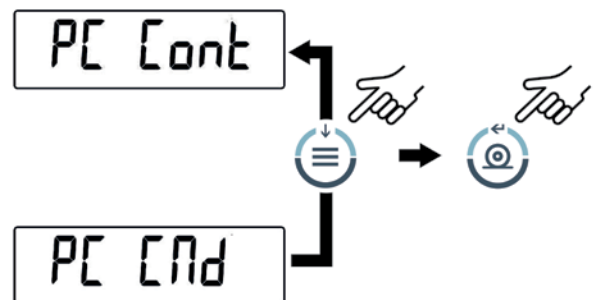Connect the balance to PC with proper cable (par. 25)

1. From zero display condition press and keep pressed the  button until the acoustic alarm gets mute, then release the button. It will be displayed the string *units*, then press the  button until it is displayed *PC-Prtr* message then confirm by pressing  button.

2. Press the  button until the *PC Cont* is displayed to choose continuous print or *PC End* to choose print to PC at user command, then press the  button to confirm the choice

3. After having selected the desired transmission mode, press the  button to go to next parameter or the  button to go to previous.

4. To escape from parameters setup menu, press the  button until the acoustic alarm gets mute, then release the button. The balance returns to normal weighing conditions.

5. The balance returns to normal weighing conditions sending the data continuously or at user command.


**NOTE:** select speed transmission (par. 12)

## 11. PRINTER SERIAL COMMUNICATION SELECTION

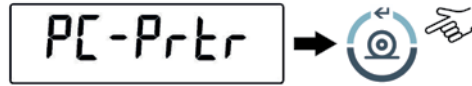Connect the balance to the printer using the proper cable (par. 25)

1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *units* will be displayed, then press  until you read the message *PC-Prtr* on display and confirm pressing  button.


2. To select the data print mode, press  until the message *Print* is displayed or *TLP50* if you want to connect the TLP50 printer for printing on labels or on paper with date and time.

3. Press  to confirm.

4. After having selected the data print mode, press  button to go to next parameter or  button to go to previous one.

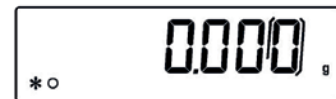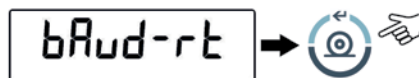
5. To escape from parameters setup menu, press  button until the acoustic alarm gets mute, then release the button. The balance returns to normal weighing conditions, ready to transmit data each time the  button is pressed.

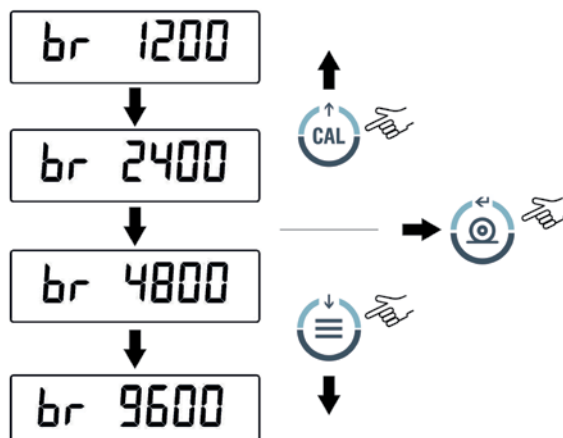
**NOTE:** elect speed transmission (par. 12).

## 12. TRANSMISSION SPEED SELECTION

1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *units* is displayed, then press  button until the message *bAud-rt* is displayed and confirm by pressing the  button.

2. Select serial data transmission speed (1200-2400-4800-9600 baud). Pressing  or  buttons it will be possible to scroll forward or backward trough the different transmission speeds, then confirm your choice by pressing  button.

3. After having selected the transmission speed you wish, press the  button to go to next parameter or  button to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm gets mute, then release the button.

5. The balance will return to standard weighing conditions.

## 13. AUTOZERO FUNCTION

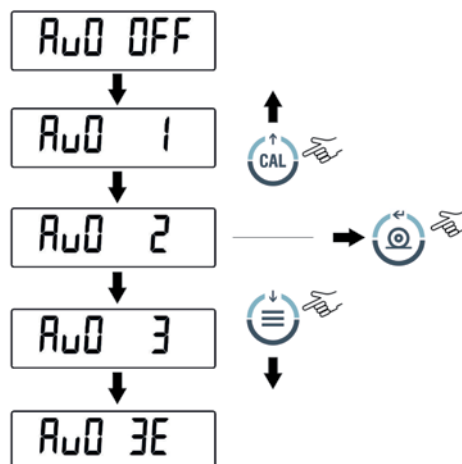
Autozero is an automatic correction of a possible zero drift.

- Au0 OFF = autozero disabled
- Au0 1 = soft autozero
- Au0 2 = medium autozero \*
- Au0 3 = heavy autozero \*
- Au0 3E = heavy autozero over all range \*      \* locked in verified balances

1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *units* is displayed, then press  button until the message *Auto 0* is displayed, then press  to confirm.

2. Pressing  or  button it will be possible to scroll forward or backward through the different autozero levels, select the one you wish and confirm it by pressing the  button.

3. After having selected the autozero level you wish, press the  button to go to next parameter or  button to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm gets mute, then release the button.


5. The balance will return to standard weighing conditions.

## 14. FILTERS SELECTION

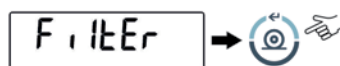
It is possible to adapt the balance to the different environment conditions thanks to the selection of three filters:

- **FILTER 1**: proportion of ingredients condition
- **FILTER 2**: stable conditions
- **FILTER 3**: unstable conditions

1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *UNIT5* is displayed, then press  button until the message *FILTEr* is displayed then confirm it by pressing the  button..

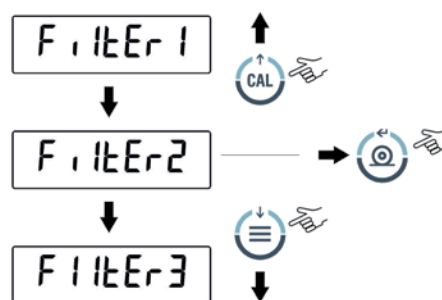
2. Pressing  or  button it will be possible to scroll forward or backward the different filtering levels, select the one you wish and then confirm it by pressing the  button.

3. After having selected the filtering level you wish, press the  button to go to next parameter or  to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm is over, then release the button.

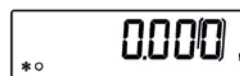5. The balance will return to standard weighing conditions.


**NOTE:** It is suggested to use FILTER 1 when proportion of ingredients must be performed

## 15. STABILITY FUNCTION

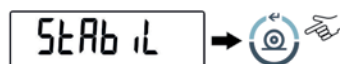
The stability symbol is displayed when the weight is stable inside a defined range

- **STAB 1** = for stable environments
- **STAB 2** = for not so stable environments
- **STAB 3** = for unstable environments

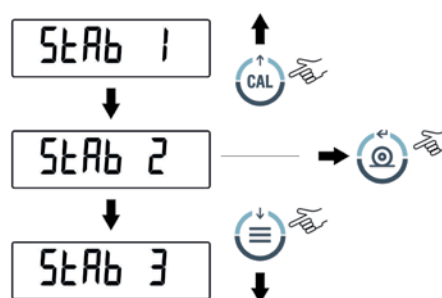1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *UNIT5* is displayed, then press  button until the message *StAbiL* is displayed, then confirm this by pressing the  button..

2. Pressing  or  button it will be possible to scroll forward or backward the different stability levels, select the one you wish and then confirm it by pressing the  button.

3. After having selected the stability level you wish, press the  button to go to next parameter or the  button to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm gets mute, then release the button.

5. The balance returns to normal weighing conditions.

## 16. DISPLAY CONTRAST REGULATION

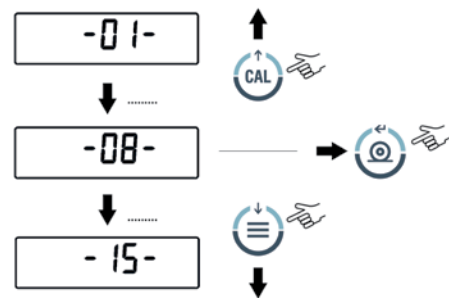
It is possible to regulate the contrast level of display to have a comfortable view of the indication at different angles of usage.

There are 15 different levels of regulation:

1. From zero condition on display, press and keep pressed the  button until the acoustic signal gets mute, then release the button. The message *units* is visualized, now press the  button repeatedly until the message *Contr* is displayed and confirm this by pressing the  button.


2. Now, pressing the buttons  or  it is possible to increment or decrement the level of the display contrast; select the one desired and confirm by pressing the  button.

3. After you have confirmed the contrast level desired, press the  button to go to the next parameter or the  button to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm is over, then release the button.

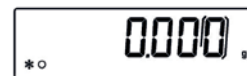5. The balance will return to standard weighing conditions.

## 17. DISPLAY BACKLIGHT SETUP

The balance display is equipped with backlight to make the indication more visible also during low light conditions.

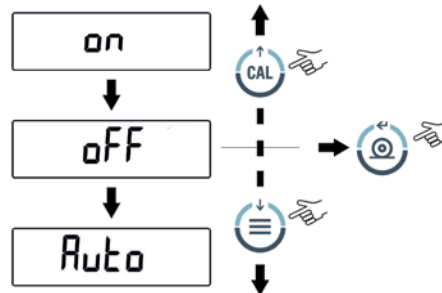
There are 3 working modes:

- **ON** = backlight always switched ON
- **OFF** = backlight always switched OFF
- **AUTO** = backlight automatically switched on during weighing operations

1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *units* is displayed, then press  button until the message *blt*, then press the  button to confirm this.


2. Pressing  or  button it will be possible to scroll forward or backward the different working modes, select the one you wish and then confirm it by pressing the .

3. After having selected the backlight working mode you wish, press the  button to go to next parameter or the  button to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm is over, then release the button.

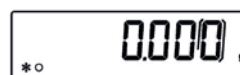5. The balance will return to standard weighing conditions.

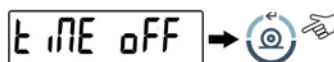
## 18. AUTO POWER-OFF FUNCTION

This function allows to activate the automatic power-off of the balance after a defined balance idle-time.

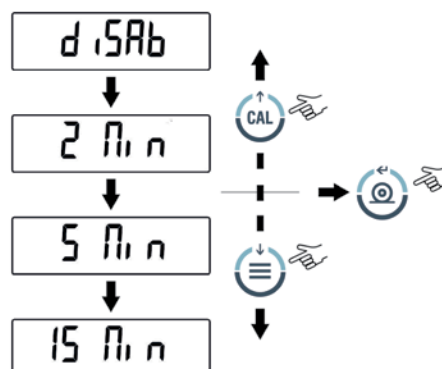
There are 4 auto power-off modes:

- **disab** = Auto power-off disabled
- **2 Min** = Auto power-off after 2 minutes of idle time
- **5 Min** = Auto power-off after 5 minutes of idle time
- **15 Min** = Auto power-off after 15 minutes of idle time

1. From zero condition on display, press and keep pressed the  button until the acoustic alarm gets mute, then release the button. The message *LINE OFF* is displayed, then press  button until the message *LINE OFF* then confirm by pressing the .

2. Pressing  or  button it will be possible to scroll forward or backward the different auto power-off modes, select the one you wish and then confirm it by pressing the .

3. After having selected the auto power-off mode you wish, press the  button to go to next parameter or the  button to go to previous one.

4. To escape from parameters setup menu, press the  button until the acoustic alarm is over, then release the button.

5. The balance will return to normal weighing conditions.

## 19. SELECTION OF WORKING MODE WITH TABLET

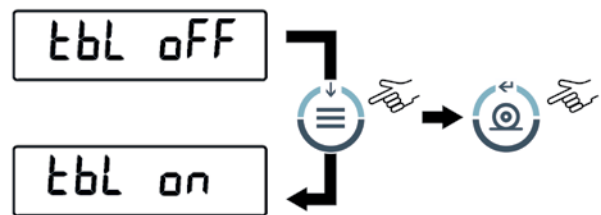
**i** NOT available for verified balance and version software RL 3.xx or R 3.xx.

Using the appropriate connection box, connect the tablet to the serial output of the balance

1. From condition of zero on display, press and hold the  button until the acoustic signal is over, then release the button. The message "unitS" is displayed, press the button  until the message *tbl Node* is displayed and confirm pressing  button

2. To choose the working mode with tablet, press the  button until you visualize the message *tbl on*, then press  to confirm

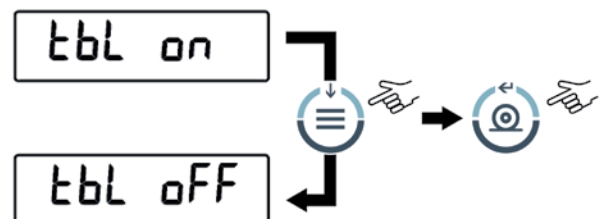
3. After you have chosen the mode "TBL ON" the balance will automatically restart and the message *TABLET* will appear on the display of balance. From this moment the interface device with the balance is the tablet.

*TABLET*

4. To escape from "TABLET" mode, press the  button, then  button.

*tbl Node*

5. Press the  button until you visualize the message *tbl off*, then press  to confirm the selection.

6. After you confirm with the  button, the balance will restart automatically and will return to standard working mode.

## 20. PIECECOUNTING FUNCTION

1. From zero condition on display, press MENU button until the message *Count* is displayed then press the  button to confirm.

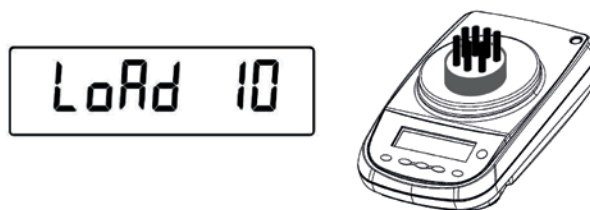
2. Select the number of pieces to put on the pan as sample by pressing  repeatedly to increase and  to decrease the number. Choice of the number of pieces (10, 25, 50, 100) depends from the weight of single pieces.

If available, load the empty container first then press  to confirm. Load on the pan the number of pieces displayed on the balance display.

Press again the  button and wait the weight to stabilize.

If the number of samples is enough (for example 10 as in figure), then this number will be displayed and it will be possible to go on by loading the pieces to count on the balance pan.

If the weight of the pieces to count is too low in comparison with balance resolution, then an error message will be displayed. In this case it is necessary to use a balance with higher resolution.

If the weight of samples is acceptable but not enough, then the message *Add* will be displayed.

Add a quantity of pieces so that to almost double the number read on display, then press the  button.

If the number of pieces is still not enough then the message *Add SNP* is displayed. Add again a quantity of pieces so that to almost double the number you read on display.


As soon as it is reached a sufficient number of pieces, this number is displayed and it will be possible to go on with count by loading the pieces count on the balance pan.

3. To escape from piececounting mode press the  button and the balance returns to normal weighing conditions.

## 20.1 VISUALIZATION OF TOTAL AND UNIT WEIGHT OF PIECES

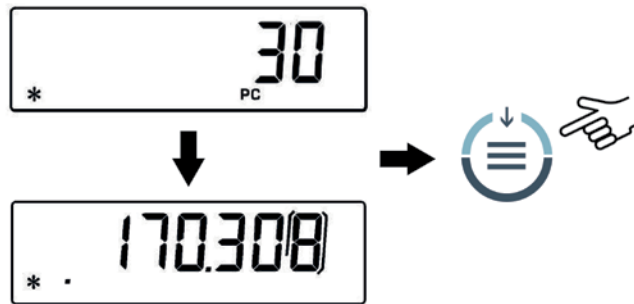1. Press  button to pass from pieces number visualization to total weight.

2. In order to display the weight of single piece from the total number of pieces press and keep pressed the  button until the acoustic alarm gets mute.

3. Press again the  button to display the number of pieces.

## 20.2 VISUALIZATION OF TOTAL AND UNIT WEIGHT OF PIECES

It is possible to insert the unit average weight of the sample, if known. In this way can be avoided the sampling of the units.

1. From zero condition on display, press the  button until the message *Count*, then press the  button to confirm. Select *NANUAL* by pressing repeatedly the  button.

2. Press the  button to confirm.

3. Insert the unit weight of a sample in grams using  and  buttons for increase and decrease the value and use the  button to skip to the others digits. For inserting the decimal point keep pressed the  button. Keeping pressed the  button is possible to delete the inserted value

4. Press the  button to confirm. If the inserted weight is 100 times lower than balance resolution, it will be displayed an error message. To escape without inserting the weight press .

5. If the weight is enough then  is displayed ; it is now possible to go on to count loading the pieces on the pan.

6. Press  button to display the total number of pieces, then press again the button to display again the number of pieces.

7. To escape from piececounting function, press  button.

## 20.3 AUTOMATIC UPDATE OF UNIT WEIGHT

After having effected the sampling, it is possible to update the average piece weight as follows:

1. Instead of loading all the pieces to count, load a number of pieces equal to the double of the ones loaded on the pan and wait for the acoustic alarm.
2. It is now possible to repeat this procedure up to a maximum of 255 pieces or to proceed to the normal counting of the pieces. This procedure allows a more accurate evaluation of the average unit weight and a better precision in the pieces count.

**NOTE:** automatic update is not active if the sampling has been effected by insertion of the average unit weight.

## 21. DENSITY DETERMINATION OF A SOLID OR A LIQUID

\* The balance must be equipped with the proper optional software (not available for all models) \*

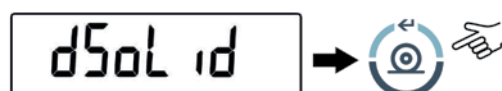
The software for density determination allows to calculate the density of a solid or a liquid through the under balance hook or through the optional hydrostatic kit.

### 21.1 DENSITY DETERMINATION OF A SOLID

1. From zero condition on display, press the  button until the message *dEnS* is displayed, then press the  button to confirm.

2. Select the function *dSol:d* through the  button and then confirm with  button.

3. The density of the liquid to use will be displayed, the default value is equal to 1.0000 (distilled water at 20°C).

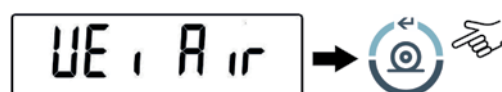
4. It is possible to set a different value using the  e  buttons for increase and decrease the value and using the  button for skip to the others digits. Keeping presses the  button is possible to delete the inserted value.

5. It is also possible to set a different value if the balance is equipped with the optional alphanumeric keyboard.

6. Once the wished value is set, press  button.

7. Now it will be asked to weigh the solid in air, confirm this with  button.

8. If necessary, effect a tare operation and load the solid. Press the  button to acquire the value. During value acquisition the message *UE, Air* will be flashing.

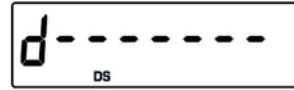
9. Now it is asked to weigh the solid inside the liquid. Effect the tare of the small basket inside liquid, immerse the solid and press . During the acquisition of the value the message *UE, LIQ* will be flashing.

10. The result of density of the solid is then displayed. Pressing the  button it is possible to print the value of density, if the balance is equipped with a printer.

11. If any error occurs, this message will be displayed:

12. Press now the  button to escape from density function, or the  button to perform another measurement.

## 21.2 DENSITY DETERMINATION OF A LIQUID

1. From zero condition on display, press the  button until the message *dEnS* is displayed, then press the  button to confirm.

2. Select the function *d Liquid* through the  button and then confirm with  button.

3. It will be displayed the value of the density of the glass float of known volume to use, the default value is equal to 3.0000.

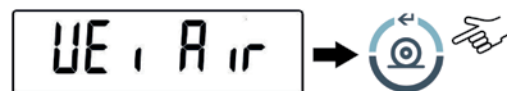
4. It is possible to set a different value using the  e  buttons for increase and decrease the value and using the  button for skip to the others digits. Keeping presses the  button is possible to delete the inserted value.

5. It is also possible to set a different value if the balance is equipped with the optional alphanumeric keyboard.


6. Once the wished value is set, press  button.

7. It is now asked to weigh the glass float in air, confirm this with  button.

8. If necessary, effect a tare operation and load the solid. Press the  button to acquire the value. During value acquisition the message *UE: Air* will be flashing.

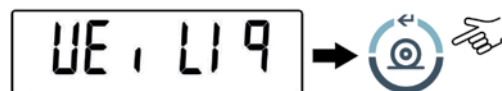9. It is now asked to weigh the glass float immersed into the liquid. Perform a tare if necessary and immerse the glass float into the liquid. Press the  button. During the acquisition of the value the message *UE: LI9* will be flashing.

10. The result of the density of the liquid is now displayed. Pressing the  button it is possible to print the value of density, if the balance is equipped with a printer.

11. If any error occurs, this message will be displayed:

12. Press the  button to escape from density function, or the  button to perform another measurement.

## 22. MAXIMUM LOAD DETERMINATION FUNCTION

**\*The balance must be equipped with the specific optional software (not available for all models)\***

The function M LOAD allows to measure the maximum load for a solid.

1. To activate the function M LOAD from zero display condition, press repeatedly the  button until the message *M LoAd* is displayed, then press the  button to confirm. When this function is activated, a tare is automatically performed.

2. Activation of this function is indicated by the visualization of the letter *M* on the last digit on the display.

3. It is now possible to determine the weight of the maximum load.

4. Then press the  button to effect another measurement.

5. When the function M LOAD is active, the calibration is disabled.

6. Press the button  to escape from M LOAD function.

## 23. PERCENTAGE WEIGHING FUNCTION

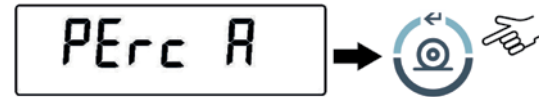
This function allows to read the weighing value as a percentage of a reference weight. The reference weight is took as 100% percentage value (factory setting). There are two modes for acquiring the reference weight: one is automatic (with the reference weight) and one manual (by inserting manually the value of the reference weight).

### 23.1 MODE WITH WEIGHT REFERENCE


1. From zero condition of the display press the  button until will be displayed the *PErc* message, then press the  button to confirm.

2. Press the  button to select the *PErc A* function and press  button to confirm.

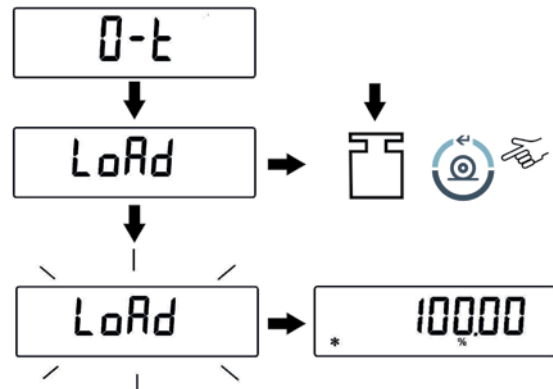3. Tare function will be performed and will be shown on the display the string *LoAd*. Then load on the weighing pan the reference and press the  button, the string *LoAd* will start to flash and once the weight is acquired will be visualized the value with the % indication.

4. Remove now the reference weight and load the sample. Is possible to read the weight in percentage now.

5. Press the  button to visualize the weight in gram and vice-versa.

6. Press the  button to escape from the percentage weight function.

**NOTE:** If the reference weight is lower than 10 digits after the acquiring sequence of the weight, the string *Error 07* will be displayed.

### 23.2 MODE WITH MANUAL INSERTION OF THE REFERENCE WEIGHT

1. From zero condition of the display press the  button until will be displayed the *PErc* message, then press the  button to confirm.

2. Press the  button to select the *PErc N* function and press  button to confirm.

3. It is possible to set a different value using the  e  buttons for increase and decrease the value and using the  button for skip to the others digits. Keeping pressed the  button is possible to delete the inserted value. The inserted value will be stored in the memory until the balance is switched off.

4. Press button when the desired value of the reference weight is inserted.

5. Load now the sample and read the percentage value.

6. Press the button to visualize the weight in gram and vice-versa.

7. Press the button to escape from the percentage weight function.

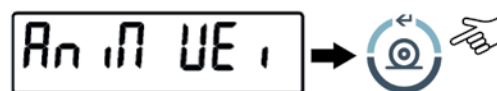
8. It is also possible to set the value if the balance is equipped with the optional alphanumeric keyboard. Press button to confirm the inserted value.

**NOTE:** If the reference weight inserted is lower than 10 digit displayed, will be shown the *Error 07*.

## 24. ANIMAL WEIGHING FUNCTION

This function allows to acquire the averaged weight of moving objects or animals for a time period that can be set.

1. From zero condition of the display press the button until will be displayed the *Anim UE* message, then press the button to confirm.

2. Set the desired period time from 5 to 90 sec. using the button to decrement and CAL to increment time. Confirm by pressing the button.

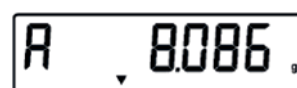
3. It is now displayed the value of current weight, effect tare if necessary, then load on the pan the sample to weight and press the button.

4. It is then displayed the count-down of the sampling time value that has been previously set.

5. When the count-down is finished, it is calculated and then displayed the value of the averaged weight measured.

6. Press once the button to start another measurement, or twice to escape this function and return to weighing mode


## 25. TOTALIZING

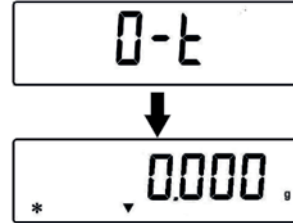This function allows to effect sum of consecutive weighings.

1. From zero condition of the display press repeatedly the  button until will be displayed the *Add* message, then press the  button to confirm.

2. The Tare is automatically effected.


3. Load the first weight to sum.

4. Confirm with the  button.

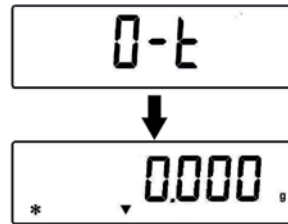5. The Tare is again effected.

6. Load the second weight to sum.

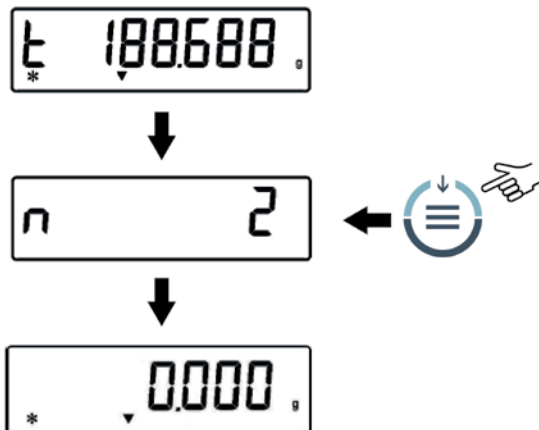
7. Confirm with the  button. Follow all the above steps for all weights to sum up to a maximum number of 99.

8. Pressing the  button it is possible to show in sequence the informations respectively of total weight, number of weighings effected and current weight.

Before the output function:

If the balance is equipped with a printer or connected to a PC, automatically after pressing the print button, will be printed or sent to PC the weigh number and its value.

After running all the desired weighs, press  to print or send to the PC the value of the total weight. The count of the weighings will then be reset and it will be possible to start over a new measure.

9. Press the  button to escape this function and return to weighing mode


## 26. THRESHOLD FUNCTION

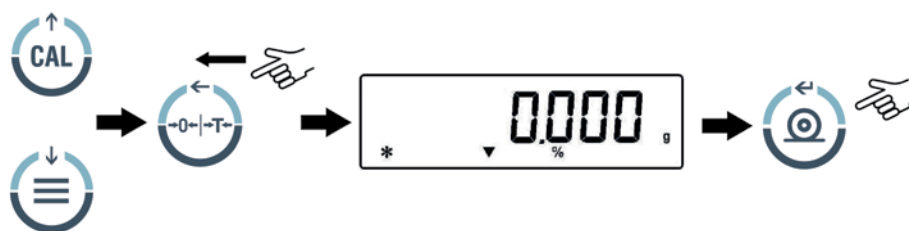Threshold function allow to determinate if the weight loaded on the pan is above or below the two threshold fixed by the user.

1. From zero condition of the display press the  button until will be displayed the string *H o L*, then press the  button to confirm.


2. Insert the lower threshold value using the  e  buttons for increase and decrease the value and using the  button for skip to the others digits. Keeping pressed the  button is possible to delete the inserted value. The inserted value will be stored in the memory until the balance is switched off.

3. Insert the higher threshold value as same as the lower value steps.

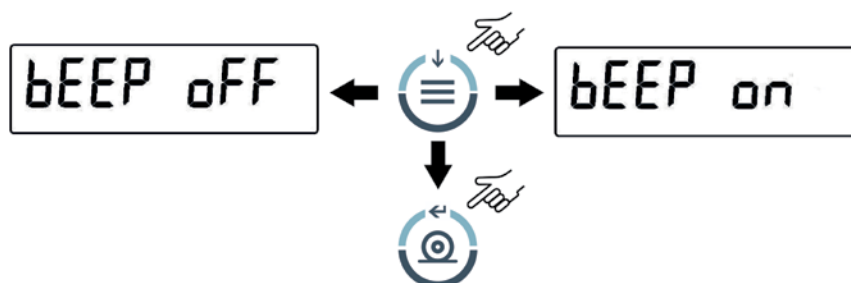4. The string *bEEP oFF* will be displayed and select by the  button if the acoustic alarm should be set or not if the weight is within the two threshold values set before. Press  button to confirm.

5. If the value of the thresholds is correctly inserted, the balance will be back to the weighing mode with the indication of the threshold state (H = higher threshold, L = lower threshold, OK= weight within the two set thresholds).

**NOTE:** If the values are not inserted correctly will be displayed the string *Error 07*.

There are 3 modes of the threshold function:

### 26.1 BOTH THRESHOLDS SET

This mode allow to set the lower and the higher threshold values and to identify the acceptable range for the weight value shown by the switching on of the symbol *OK* and by the acoustic alarm is activated. When the value of the weight is under the lower threshold value set will be displayed the symbol *L* while if the weight is over the higher threshold value set will be displayed the symbol *H*.

### 26.2 LOWER THRESHOLD SET ONLY

Setting only the lower threshold and setting to zero the higher threshold, will be displayed the symbol *OK* and eventually the acoustic alarm if activated whenever the loaded weight is higher than the lower threshold value set. When the value of the weight is under the lower threshold set the symbol *L* will be displayed.

## 26.3 HIGHER THRESHOLD SET ONLY

Setting only the higher threshold and setting to zero the lower threshold, will be displayed the symbol  $OK$  and eventually the acoustic alarm if activated whenever the loaded weight is lower than the higher threshold value set. When the value of the weight is above the higher threshold set the symbol  $H$  will be displayed.

# 27. RS232 INTERFACE FEATURES

## 1. General features

The balance transmits the value visualized on the display following serial RS232C standard, allowing to print the value of weight to a PC monitor or to a serial printer. In the case of connection to a PC, it will be possible to select the transmission in continuous mode or transmission at user command through pressing of the  button (as described at par.10). The balance is also capable of receiving commands,

always through the standard RS232C, that allow performing all the functions available through the keyboard of PC itself. The speed of transmission and reception can be selected, as described previously (see par.12), to 1200, 2400, 4800, e 9600 baud. The character format is of 8 bit preceded by one bit of start and followed by a bit of stop. Parity is not considered.

## 2. Selection of interface for PC

Selecting the transmission to PC (personal computer IBM compatible), it is achieved a continuous transmission output, at the same rate of weight update on display of the balance. It is possible to perform all the functions of the balances directly from the computer keyboard, transmitting to the balance the ASCII codes as in the table below. The connector to use for connection to PC is the number 2 with reference to fig.2 of paragraph 5 of this manual.

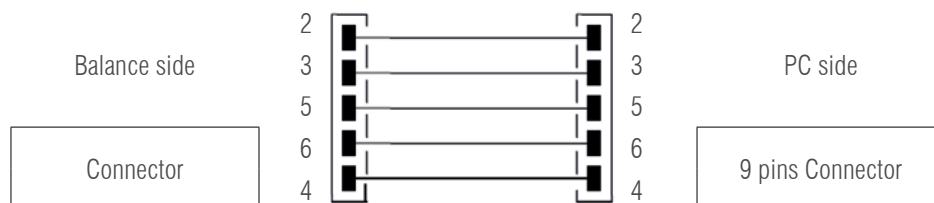
CODE		1st FUNCTION (SINGLE PRESS)
ASCII	HEX	
T	54	TARE
C	43	CALIBRATION
E	45	ENTER
M	4D	MENU
O	4F	ON/OFF

CODE		2nd FUNCTION (PROLONGED PRESS)
ASCII	HEX	
t	74	TARE
c	63	CALIBRATION
e	65	ENTER
m	6D	MENU
o	6F	ON/OFF

Selecting the transmission to PC at user command, it is achieved a transmission output only when the  button is pressed, also in this case it is possible to perform all the functions of the balance directly from the keyboard of the computer, sending to the balance the ASCII codes in the table above.

## 3. Connection of the balance to PC

To receive/transmit data, link the connector (n.2 of fig.2 at par.5) of the balance to the serial port of your Personal Computer as shown below:

#### 4. Transmission format

String transmitted is composed by the following 14 characters:

- First character: weight sign (blank or -)
- second/ninth character: weight or other data
- tenth/twelfth character: weight unit symbol
- thirteenth character: stability indicator
- fourteenth character: carriage return
- fifteenth character: line feed

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Eventual non-significative zero are put as spaces.

In the following tables the various transmission formats are shown:

WEIGHING MODE (VALID BOTH FOR CONTINUOUS TRANSMISSION AND TRANSMISSION AT USER COMMAND)

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15
Sign	Weight							Weight unit			Stability	CR	LF	

DENSITY MODE (ONLY IN TRANSMISSION AT USER COMMAND MODE)

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17
d	=	Density value						Space	Weight unit				CR	LF		

PIECECOUNTING MODE (ONLY IN TRANSMISSION AT USER COMMAND MODE)

- Number of pieces

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	
Pcs			:	spaces				Number of pieces								

- Total weight of pieces

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18	19	20
Weight						:	spaces	Value of weight							space	g	space	S	

- Average unit weight of pieces

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
PMU			:	spaces				Value of weight							space	g	

PMU stands for average unit weight

PERCENTAGE WEIGHT MODE (ONLY IN TRANSMISSION AT USER COMMAND MODE)

- Percentage

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Perc			.	space				Percentage							space	%	

- Weight

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Weight				space				Weight value							space	%	

ANIMAL WEIGHING MODE (ONLY IN TRANSMISSION AT USER COMMAND MODE)

• Time

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Time				space	=	space			Time value			Sec		space			

• Averaged Weight

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Ave		.	=	space				Averaged weight value					space		g		

TOTALIZING (ONLY IN TRANSMISSION AT USER COMMAND MODE)

• Weighing

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Weighing number		.	space			Value of weight						space		g			

• Weighs total

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
S	space	=	space			Value of weight						space		g			

THRESHOLD FUNCTION (ONLY IN TRANSMISSION AT USER COMMAND MODE)

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18	19	20
Weight						:	minus sing if negative	Value of weight						space		g			

If Low

1	2	3	4	5
- Low -				

Else Hight

1	2	3	4	5
- Low -				

Else ok

1	2	3	4	5
- Ok -				

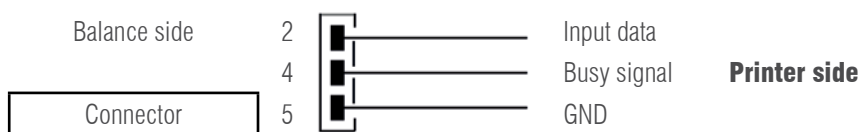
**5. Selection of interface to printer**

Selecting the PRINTER mode, the serial output of the balance is set to work with serial printers.

In this case the printing is effected only when the  button is pressed and with stable weight. If the stability is not reached within ten seconds, the message *ERR05* is displayed preceded by a short acoustic signal and the value of the weight is not sent to the printer.

**6. Connection of the balance with the serial printer**

Connect a serial printer to the connector (n.2 of fig.2 at par.5) to the balance as shown in the following scheme:

If the optional TLP50 printer is used, it will be possible to print both in continuous-mode and in labels-mode with the following formats:

Weighing mode and maximum load mode

12-02-2008	12:00
Weight:	22.000 g

Piececounting mode

12-02-2008	12:00
Pcs	100
Weight:	300.000g
PMU:	3.000g

Density determination mode

12-02-2008	12:00
d=	2.80066 g/cm <sup>3</sup> d

Percentage weight mode

12-02-2009	12:00
Perc.	100.0%
Weight:	300.000 g

Animal weighing mode

12-02-2009	12:00
Time =	6 Sec
Ave. =	59.446 g

Totalizing (sum of weights) mode

12-02-2010	12:00
1.	16.589 g
2.	17.226 g
...	
99.	
.....	
S=	33.815 g

Threshold function

12-02-2013	12:00
Weight:	0.00g
	-LOW-

12-02-2012	12:00
Weight:	49.20g
	- OK -

12-02-2011	12:00
Weight :	249.42g
	-HIGH-

**7. Map of RS232 interface Connector (n.2 fig.2 par.5)**

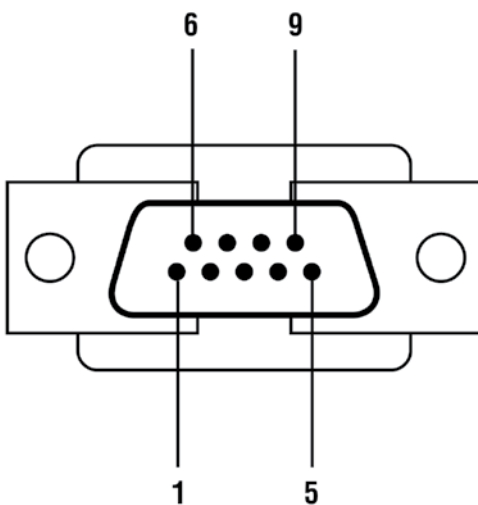
Fig. 1

CONNECTOR 1: CONNECTION FOR KEYPAD, PC AND PRINTER

- pin 1 = Power +5V for keyboard
- pin 2 = Tx signal
- pin 3 = Rx signal
- pin 4 = busy signal
- pin 5 = Gnd
- pin 4-6 = connected to each other for connection to PC

Fig. 2

## 28. ERROR CODES

ERROR MESSAGE ON DISPLAY	MEANING	POSSIBLE SOLUTIONS
<i>Erro1</i>	Weight not stable after operation of tare	Protect the balance from air flows or from vibrations of the working table
<i>Erro2</i>	impossible to start the calibration due to instability of the balance	Protect the balance from air flows or from vibrations of the working table.
<i>Erro3</i>	calibration weight not correct or balance unstable	Calibrate with correct weight or protect the balance from environment disturbs
<i>Erro4</i>	weight of samples for the piececounting function not adequate or unstable	Select a bigger number of samples or protect the balance from vibrations.
<i>Erro5</i>	impossible to print / weight unstable	Protect the balance from environment disturbs
<i>Erro6</i>	Weight cannot get stable in density mode	Protect the balance from environment disturbs
<i>Erro7</i>	error in inserting the data	
<i>Erro8</i>	Anomaly on autocalibration motor	Contact service staff
<i>unLoAd</i>	weight loaded on the pan or pan not positioned properly	Remove the weight from the pan or position properly the pan and underpan.
<i>CAL but</i>	the balance requires to be recalibrated	Unload weights, if any, on the pan, and press the CAL button
	Over-range condition	Unload the weights loaded on the pan
	Under-range condition	Place properly pan and underpan

## 29. MAINTENANCE AND CARE

Regular maintenance of yours balance guarantee accurate measurements.

### 1) Cleaning

Before cleaning the balance unplug the power supply of the balance from the voltage supply of your room. Do not use aggressive cleaning product (as solvents or similar), use a humid towel with soft detergent, Avoid liquids to go inside the instrument during the cleaning. Wipe the balance with a soft towel. Parts of samples or powder can be removed using a brush or vacuum cleaner.

### 2) Safety Checks

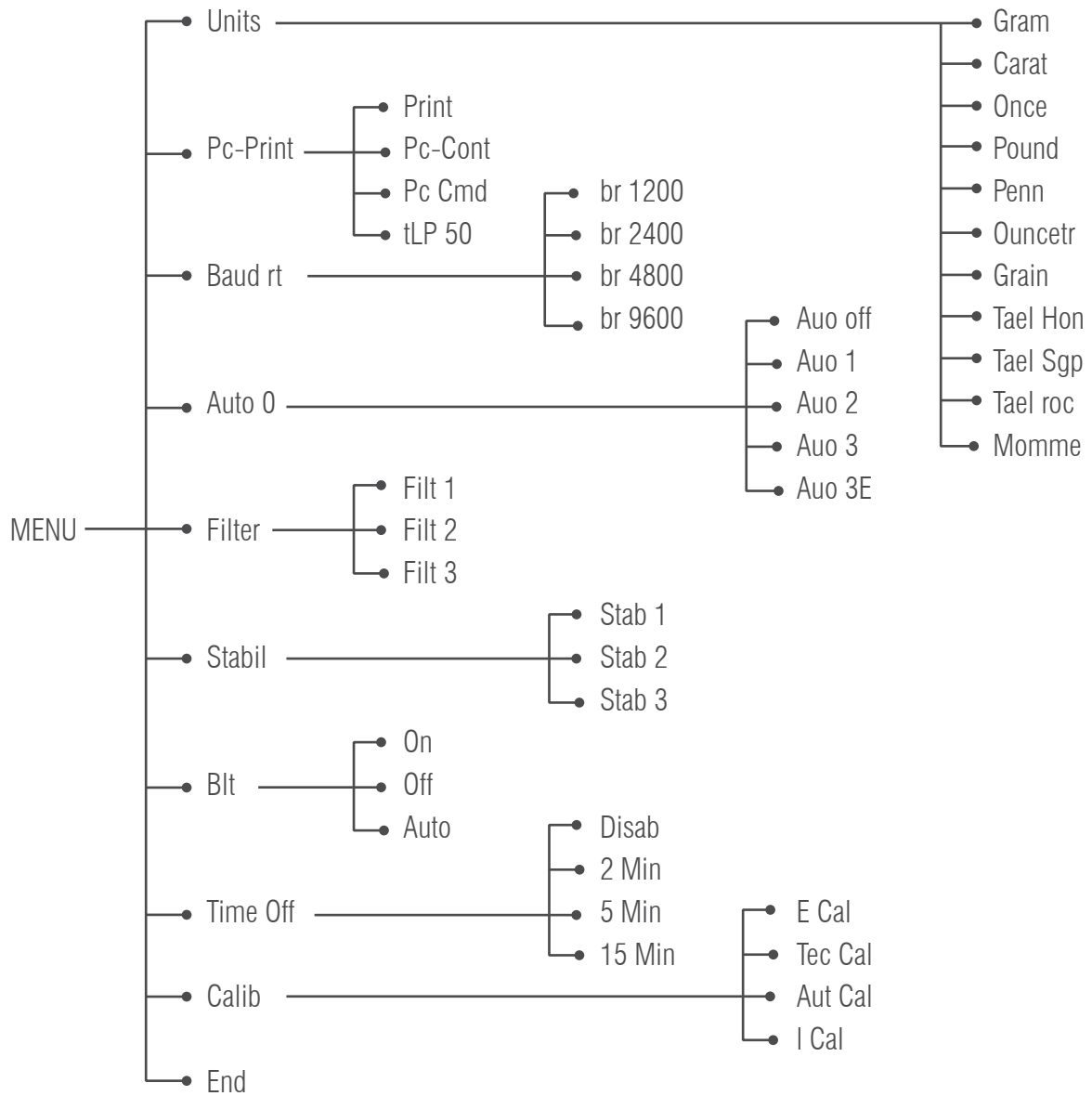
Safety of the instrument is no more guaranteed when:

- balance power supply is clearly damaged
- balance power supply is not working anymore
- balance power supply is stored for long time in hard environment conditions.

In these instances refer to the assistance centre where specialized technician will make reparations to bring back the instrument in the safety conditions eventually.


### 30. QUICK GUIDE TO BALANCE PARAMETERS SETUP

- To enter the balance parameters setup menu, press and keep pressed the  button until the acoustic alarm gets mute.
- Use then the  button to go to next parameter, use the  button to go to previous and the  button to confirm the choice.
- To escape from menu, press and keep pressed the  button until the acoustic alarm gets mute.

## 31. QUICK GUIDE TO THE USE OF THE BALANCE PROGRAMS

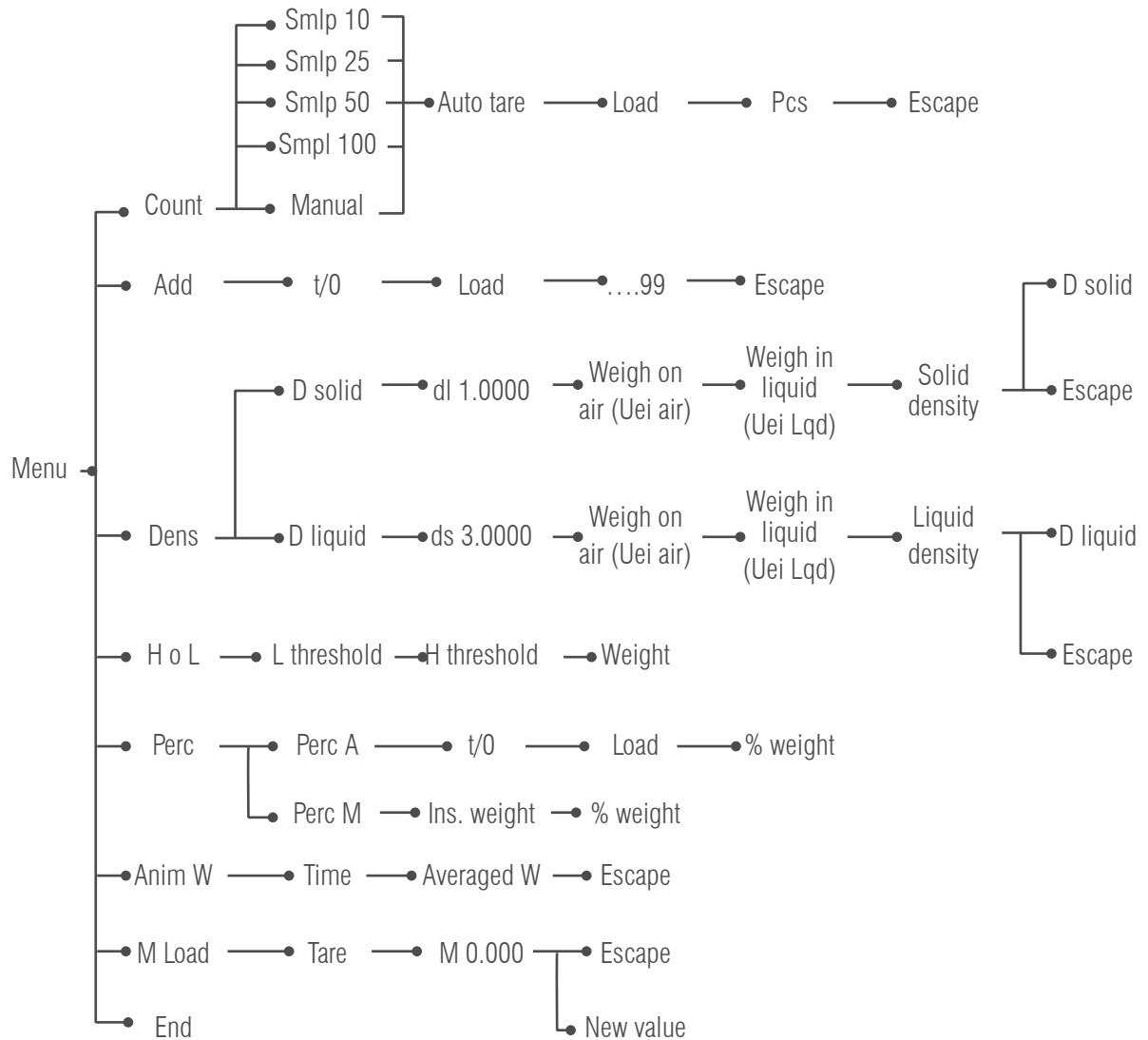- To enter the balance programs menu, press the  button.
- Use then the  button to go to next parameter, use the  button to go to previous and the  button to confirm the choice.
- To escape from menu, press and keep pressed the  button until the acoustic alarm gets mute.

## 32. BALANCE TECHNICAL CHARACTERISTICS

The models below indicated are all for internal use. Max altitude use: 4000m; Pollution grade: 2; Overvoltage category: II

Power supply included:	INPUT: Switching 100-240Vac~ 50/ 60Hz, OUTPUT: 24V DC 550mA, Max absorbed power 13.2VA
Environment condition adaption:	Selectable filters
Autozero:	Selectable from Menu
Serial output:	RS232C
Operating temperature range:	+5°C - +35°C

Power supply included:	INPUT: Switching 100-240Vac~ 50/ 60Hz, OUTPUT: 9V DC 1000mA, Max absorbed power 9VA
Environment condition adaption:	Selectable filters
Autozero:	Selectable from Menu
Serial output:	RS232C
Operating temperature range:	+5°C - +35°C

## 33. WARRANTY

We guarantee that under proper using situation. We provide one year repairing service include material and technical support after selling date. In Guarantee period, if machine broken or damage because of material or techniques. We will repair or replace the problem parts which has been proved. Please contact our Local office when machine need repairing.

The Guarantee Card will be inefficacy with wrong operating and not according as the operating manual. The Guarantee Card will be inefficacy with any damage or broken by unauthorized person's repairing or replacement.

We are not in charge with apparent or intentional disobeying the guarantee rule which cause machine any relevant or accidently broken.

## 34. EQUIPMENT DISPOSAL

This equipment is marked with the crossed out wheeled bin symbol to indicate that this equipment must not be disposed of with unsorted waste. Instead it's your responsibility to correctly dispose of your equipment at lifecycle -end by handling it over to an authorized facility for separate collection and recycling. It's also your responsibility to decontaminate the equipment in case of biological, chemical and/or radiological contamination, so as to protect from health hazards the persons involved in the disposal and recycling of the equipment.

For more information about where you can drop off your waste of equipment, please contact your local dealer from whom you originally purchased this equipment.

By doing so, you will help to conserve natural and environmental resources and you will ensure that your equipment is recycled in a manner that protects human health.

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marca propiedad de | est une marque de | trade mark propriety of: